



Laboratory Equipment Manufacturer
www.mrclab.com



INSTRUCTION MANUAL FOR

Circulated Water Bath

WBL-Series




PLEASE READ THIS MANUAL CAREFULLY BEFORE OPERATION

3, Hagavish st. Israel 58817 Tel: 972 3 5595252, Fax: 972 3 5594529 mrc@mrclab.com

MRC.VER.02-10.10

Name and Function of Control panel

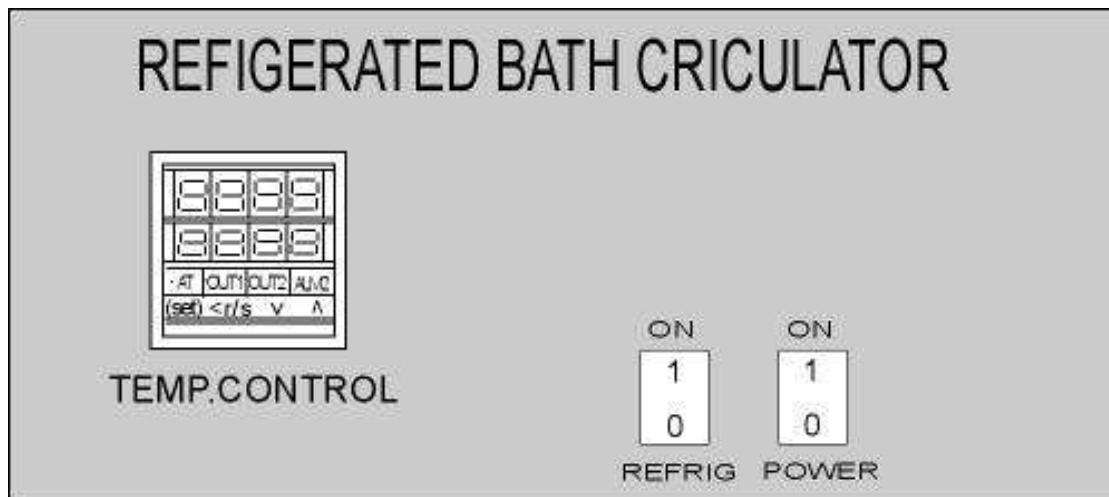

1. **Power supply switch.**

Plug in the power supply switch to its proper Outlet.

Push power control to ON position.

2. **Temperature control.**


Set the temperature to the required temperature rate.

Press the set key to enter SV mode, the digit which lights brightly is settable.

3. **Refrigerator control**

starts to cool-down minutes after turning on.

Turn on below the room temperature, avoid turning to over 40 °C room temperature as this may reduce the life span of the compressor.

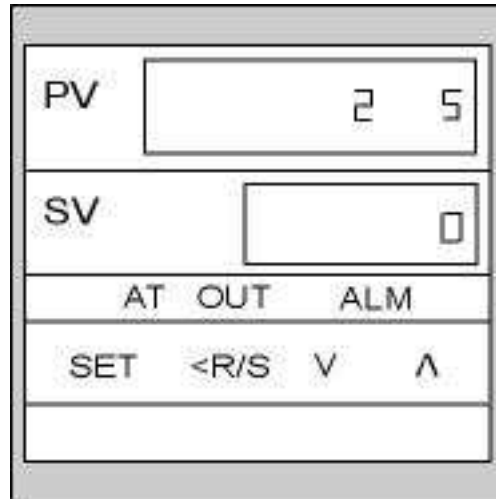

PV- display of actual temperature inside cabinet.

SV- Set-up temperature indicator.

AT- Automatic Calculation P.I.D

OUT-Control system output indicator

ALM- Over-temperature protection indicator. Once the indicator is lit, it means the temperature is over the set-up temperature range for protection.

SET-If there is a temperature change, push this button first. Then there is a 4-digit number shown on SV, and one of the lamps will be lit. The lamp can be used to set up the temperature. To push ◀ button to move left, and ▲ button to increase, ▼ button to decrease.

R/S- When using left arrow button to set up temperature, one of the lamp will be lit, and you could use this lamp to shift the position.

▼ - button to decrease set-up temperature: to decrease set-up temperature, press this button. Once completed, press SET button again.

▲ - button to increase set-up temperature: to increase set-up temperature, press this button. Once completed, press SET button again.